

www.precellys.com

NEW GENERATION Tissue homogenizer

SAMPLE PREPARATION & CELL LYSIS



precellys 24
lysis & homogenization


Transport Box

The Precellys 24 box has been especially designed to protect the equipment during transport. Whether it is by car, train, or even by plane the new transport box will protect the machine that demands a strong box that can protect its internal engine and delicate parts. With its new design the transport box can protect the Precellys from even the toughest transports. The box has four locking points, four detachable wheels, and a depressurization valve for airplane transportation.

- **Shock Resistant**
- **Easy to Use/ Transport**
- **Water Resistant and Secure**

➤ SPECIFICATIONS

- Dimensions: 52.1 x 52.1 x 48.9 cm
- Reference: 03119.870.NC005
- Weight: 13.2 kg

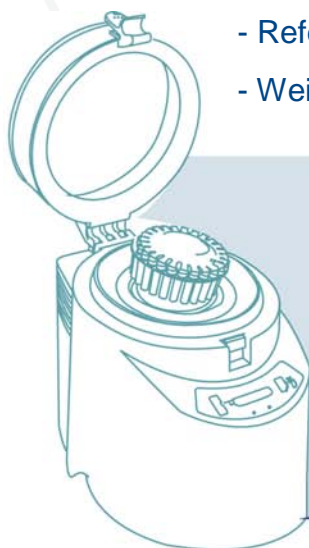

www.precellys.com
precellys@bertin.fr

Phone : +33(0)139 306 118
Fax : +33(0)139 306 185

bertin
TECHNOLOGIES
mastering innovation

For more details, please contact
precellys@bertin.fr or visit our website



www.precellys.com

NEW GENERATION Tissue homogenizer

SAMPLE PREPARATION & CELL LYSIS




precellys 24
lysis & homogenization

Transport Box Instructions

STEP #1

If needed the wheels can be found in the storage in the upper right corner of the box. To attach the wheels turn the box upside-down (with out the machine inside) and slide the wheels in their appropriate spots.

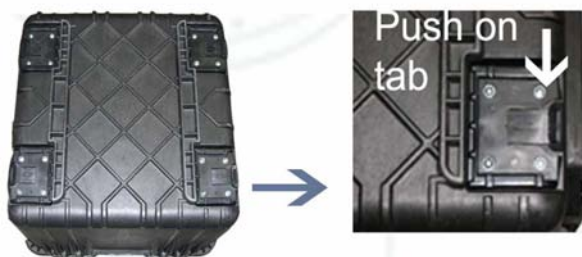

Make sure if transporting by plane, train, or automobile to remove all wheels. To remove wheels push down on the tabs and slide wheels away from box.

STEP #2

In order to make sure the machine is secure inside of the transport box, remove the white tube block plate. In addition place the protective blue foam provided over the tube receptacle. Close the hood of the machine, and remove the power cord. For added security please secure the hood with tape to ensure it does not come unfastened.



STEP #3

Unlatch the six latches found on the front, right, and left side



Place the machine facing away from you in the space provided. Make sure the machine is properly resting so that the foot rests are suspended in air.

STEP #4

Refasten all six latches



The machine is now ready for transportation.

www.precellys.com
precellys@bertin.fr

Phone : +33(0)139 306 118
Fax : +33(0)139 306 185

bertin
TECHNOLOGIES
mastering innovation

For more details, please contact
precellys@bertin.fr or visit our website

